AccuPaste™

Silver Conductive Paste

Cat. No. TC-4000-0~3

AccuPaste™ Silver Conductive Paste

User Guide

TC-4000-0~3

Version No.: 0 (2022-07-01)

Please read all the information in booklet before using the unit

BIONEER Corporation

**8-11, Munpyeongseo-ro, Daedeok-gu, Daejeon
34302, Republic of Korea**

Tel: 1588-9788

Email: sales@bioneer.co.kr

www.bioneer.com

Intended Use

AccuPaste[™] Silver Conductive Paste is developed and supplied for research purposes only. Certain applications possible with this kit may require special approval by appropriate local and/or national regulatory authorities in the country of use.

Safety Warning and Precaution

Wear appropriate protection when handling any irritant or harmful reagents. The use of a laboratory coat, protective gloves and safety goggles are highly recommended. For more information, please consult the appropriate Material Safety Data Sheet (MSDS).

Warranty and Liability

All BIONEER products undergo extensive Quality Control testing and validation. BIONEER guarantees quality during the warranty period as specified, when following the appropriate protocol as supplied with the product. It is the responsibility of the purchaser to determine the suitability of the product for its particular use. Liability is conditional upon the customer providing full details of the problem to BIONEER within 30 days.

Quality Management System ISO 9001 Certified

Every aspect of our quality management system from product development, production to quality assurance and supplier qualification meets the world-class standards.

Trademark

AccuPaste[™] is a trademark of BIONEER Corporation.

Copyright

Copyright 2022 BIONEER Corporation. All Rights Reserved.

Notice

BIONEER Corporation reserves the right to make corrections, modifications, improvements and other changes to its products, services, specifications or product descriptions at any time without notice.

Contents

Product Information	1
Components	1
Storage	1
Specifications	1
Introduction	2
Product Description	2
Features & Benefits	3
Experimental Procedures	4
Ordering Information	7
Explanation of Symbols	7

Product Information

Components

Components	Amount			
	TC-4000-0	TC-4000-1	TC-4000-2	TC-4000-3
AccuPaste™ Silver Conductive Paste	50 g	100 g	500 g	1 kg

* **Note:** Please inquiry for bulk orders.

Storage

AccuPaste™ Silver Conductive Paste must be stored in a sealed condition and at room temperature. If the storage temperature drops below zero degree, the product can become frozen. If the storage temperature rises over 90°C the composition ratio of paste can change due to the evaporation of solvent.

Specifications

Division	Standard	Note
Components	Ag powder	-
Exterior	Silver	Visual Measurement
Operating Temperature (°C)	-20 ~ 300°C	-
Viscosity(cP)	200,000	Rotary rheometer
Drying Conditions	30 minutes at 300°C	Hot Air Dry
Specific resistance ($\mu\Omega\cdot\text{cm}$)	20	ASTM D991
Storage Conditions	At room temperature	Sealed Storage

Introduction

Product Description

AccuPaste™ Silver Conductive Paste is composed of high performance silver flake powder. A coating layer maintains heat-resistance property under high temperature environment (maximum of 300°C) and very low electrical volume resistivity, unlike general conductive paste. Application of AccuPaste™ Silver Conductive Paste can be divided into printed electronics and as ESD/EMI shielding material due to its high electrical conductivity. Therefore, AccuPaste™ is applicable to manufacture premium quality products requiring operation of high volume resistivity under high temperature condition, with a slim design. Recently, there are difficulty implementing high performance conductive paste with heat resistance due to limitation of the materials and thus the demand for high-performance new materials has increased. AccuPaste™ satisfies these requirements, and the demand.

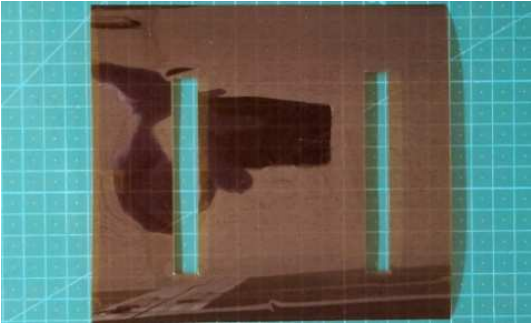
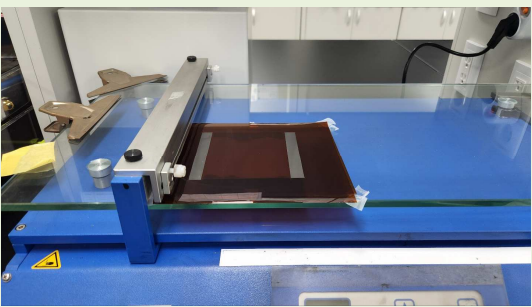
Figure 1. AccuPaste™ Silver Conductive Paste.


Features & Benefits

- Great performance
Conductive paste shows high performance using silver powder as its component.
- High thermal stability
Conduction characteristics is maintained without separation of coating film even at the maximum temperature of 300°C.
- Comprehensive compatibility
Shows good adhesion to various materials like PI(Poly Imide) film, ceramic glass, metallic substrate, etc.
- Low sheet resistance
Specific resistance is about 20 $\mu\Omega\cdot\text{cm}$.
- High electric conductivity
Applicable as electrode material with its good electric conductivity.


Experimental Procedures

The table below shows a method to use *AccuPaste™* as a pair of electrodes using the process of printed electronics. The method of its usage is not limited to the following example, and can be appropriately applied to other fields and processes.

Steps	Procedure Details
	<ol style="list-style-type: none"> 1. Prepare a sample surface to apply <i>AccuPaste™</i> to form a pair of electrodes.
	<ol style="list-style-type: none"> 2. Prepare a stencil to form electrodes with <i>AccuPaste™</i> Silver Conductive Paste by printed electronics method.
	<ol style="list-style-type: none"> 3. Place the stencil on top of the sample, parallel to the coating direction of the coating equipment.

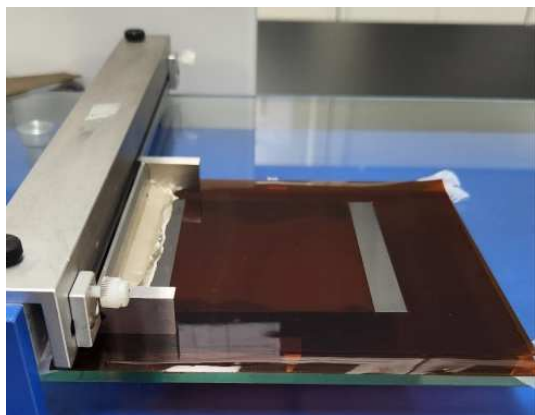4. Place a doctor blade on the sample and mount adequate amount of *AccuPaste™* Silver Conductive Paste on the doctor blade.

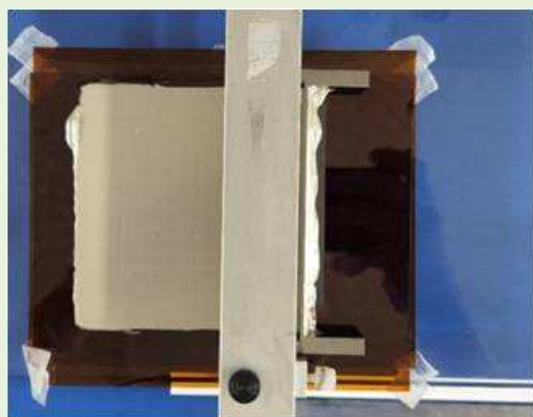5. Start to coat.

6. Uncover the stencil from the sample.

7. Cure the sample in 300°C.

Ordering Information

Description		Cat. No
AccuPaste™ Silver Conductive Paste	50 g	TC-4000-0
	100 g	TC-4000-1
	500 g	TC-4000-2
	1 kg	TC-4000-3

* **Note:** For a more detailed information of our products or to place an order, please email nano-support@bioneer.com.

Explanation of Symbols

 Batch Code	 Catalog Number	 Caution	 Consult Instructions For Use
 Manufacturer	 Research Use Only		

AccuPaste™ Silver Conductive Paste

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BIONEER Corporation - HQ
BIONEER Inc. - USA Branch
BIONEER Corp. - European Branch